
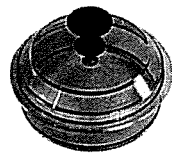


HG-4L*

Stainless Steel Rotor Body
 Aluminum Swinging Buckets
 4-Place, 1000 ml
 Tube Dimensions: 98 x 175 mm
 Maximum Speed: 5250 rpm
 Rotor Mass: 18.4 kg

| | |
|--------------|----------|
| Max. RCF: | 7120 g |
| Max. Radius: | 23.11 cm |

*Requires Conversion Kit 12300 for use in RC-3B, RC-3C.

| | |
|--|--|
| <p>Catalog No. 44000</p> <p>Includes: HG-4L Rotor Rotor Buckets, Anodized Aluminum, (2pkgs, 2/pkg) 44040 Rotor Cleaning Kit 12259 Balancing Discs, Rubber (1 pkg, 6/pkg) 00335 Instruction Manual 44039</p> <div style="text-align: center;">  <p>00335</p> </div> | <p>Accessories</p> <p>Rotor Extractor 01052 Conversion Kit For use in RC-3B and RC-3C Centrifuges 12300 Bucket Cover, Polysulfone 00395, (For 1 ml to 50 ml Tubes ONLY. CANNOT BE USED WITH STACK ADAPTER 00858) Adapter Removal Post 00392 Wire Carrying Handle for Adapters 00390, 00384, 00385. 00680</p> <div style="text-align: center;">  <p>00395</p> </div> |
|--|--|

TUBES, BOTTLES FOR HG-4L

| CAPACITY (ml) | | DESCRIPTION | MAX SPEED (rpm) | DIMENSIONS (mm) | TUBE CAT. NO. (Qty/Pkg) | COVER ASSEMBLY | | ADAPTER REQ'D | | NOTES |
|---------------|------|---|-----------------|-----------------|-------------------------|----------------------|---------------------------|--------------------------------------|---------------|---|
| NOMINAL | FILL | | | | | CAT. NO. (Qty/Pkg) | DESCRIPTION | CAT. NO. (Qty/Pkg) | NO. OF PLACES | |
| 1000 | 1000 | Polycarbonate, Oak Ridge Bottle | 5250 | 98 x 175 | 03947 (4) | 03280 (12) | Polypropylene, Sealing | — | — | |
| 1000 | 1000 | Polypropylene, Bottle | 5250 | 98 x 175 | 03695 (4) | Included with Bottle | Polypropylene, Nonsealing | — | — | |
| 450 to 600 | — | Blood Bag, Triple, Quad, and 600 ml Pack. | — | — | — | — | — | 00512 (1) | 1 | |
| 450 | — | Blood Bag, Single and Double | — | — | — | — | — | 00511 (1) | 1 | |
| 500 | 425 | Polycarbonate, Oak Ridge Bottle | 5250 | 70 x 160 | 03944 (6) | 03280 (12) | Polypropylene, Sealing | 00444 (1) 00511 (1) | 1 1 | Adapters 00444 and 00511 are both required. Nonsealing Cap 03928 (12/pkg) available. |
| 500 | 425 | Polypropylene, Oak Ridge Bottle | 5250 | 70 x 160 | 03943 (6) | 03280 (12) | Polypropylene, Sealing | 00444 (1) 00511 (1) | 1 1 | Adapters 00444 and 00511 are both required. Nonsealing Cap 03928 (12/pkg) available. |
| 315 | 360 | Stainless Steel Bottle | 5250 | 61 x 153 | 00522 (1) | 00518 (1) | Stainless Steel, Sealing | 00443 (1) 00511 (1) | 1 1 | Adapters 00443 and 00511 are both required. Requires Wrench 01014 to seal bottle. |
| 285 | 330 | Stainless Steel Bottle | 5250 | 61 x 140 | 00530 (1) | Included with Bottle | Stainless Steel, Sealing | 00443 (1) 00511 (1) | 1 1 | Adapters 00443 and 00511 are both required. |

(continued on next page)

TUBES, BOTTLES FOR HG-4L (cont.)

| CAPACITY (ml) | | DESCRIPTION | MAX SPEED (rpm) | DIMENSIONS (mm) | TUBE CAT. NO. (Qty/Pkg) | COVER ASSEMBLY | | ADAPTER REQ'D | | NOTES |
|---------------|--|---------------------------------|-----------------|-----------------|-------------------------|----------------------|---------------------------|-------------------------------------|---------------|---|
| NOMINAL | FILL | | | | | CAT. NO. (Qty/Pkg) | DESCRIPTION | CAT. NO. (Qty/Pkg) | NO. OF PLACES | |
| 250 | 280 | Polysulfone, Oak Ridge Bottle | 5250 | 61 x 124 | 03941 (12) | 03278 (12) | Polypropylene, Sealing | 00443 (1) 00511 (1) | 1 1 | Adapters 00443 and 00511 are both required. Nonsealing Cap 03927 (12/pkg) available. |
| 250 | 280 | Polypropylene, Oak Ridge Bottle | 5250 | 61 x 124 | 03937 (12) | 03278 (12) | Polypropylene, Sealing | 00443 (1) 00511 (1) | 1 1 | Adapters 00443 and 00511 are both required. Nonsealing Cap 03927 (12/pkg) available. |
| 250 | 280 | Polycarbonate, Oak Ridge Bottle | 5250 | 61 x 124 | 03939 (12) | 03278 (12) | Polypropylene, Sealing | 00443 (1) 00511 (1) | 1 1 | Adapters 00443 and 00511 are both required. Nonsealing Cap 03927 (12/pkg) available. |
| 250 | — | Conical Tissue Culture Bottle | — | 60 x 177 | — | — | — | 00443 (1) 00511 (1) | 1 1 | Adapters 00443 and 00511 are both required. Corning adapter 25351-C must be purchased from Corning Glass. |
| 200 | 200 | Nunc Tissue Culture Bottle | 5200 | 60 x 135 | 03094 (4) | — | — | 00443 (1) 00511 (1) 03095 (1) | 1 1 1 | All three adapters are required. |
| 200 | 350 | Stainless Steel Tube | 5250 | 61 x 124 | 00521 (1) | — | — | 00443 (1) 00511 (1) | 1 1 | Adapters 00443 and 00511 are both required. |
| 150 | 179 | Corex® Bottle | 5000 | 53 x 132 | 03229 (1) | Included with Bottle | Polypropylene, Nonsealing | 00372 (1) 00443 (1) 00511 (1) | 1 1 1 | One each of Adapters 00372, 00443, and 00511 are required per bottle. |
| 125 | 150 | Corex® Bottle | 5000 | 53 x 118 | 03231 (1) | Included with Bottle | Polypropylene, Nonsealing | 00372 (1) 00443 (1) 00511 (1) | 1 1 1 | One each of Adapters 00372, 00443 and 00511 are required per bottle. |
| 100 | 134 | Pyrex® Bottle | 3000 | 44 x 137 | 03230 (1) | Included with Bottle | Polypropylene, Nonsealing | 00371 (1) 00443 (1) 00511 (1) | 1 1 1 | One each of Adapters 00371, 00443, and 00511 are required per bottle. |
| 50 | Accepts all Tubes, Bottles, and Adapters for SS-34 and SA-600 Rotors; requires Adapter 00894 (1/pkg, 5-place) PLUS any adapter specified in SS-34/SA-600 tube chart. | | | | | | | | | |
| 50 | — | Tissue Culture Tube | — | 30 x 115 | — | — | — | 00390 (1) | 4 | |
| 16 | Accepts all Tubes, Bottles, and Adapters for SM-24 Rotor; requires Adapter (00892) (1/pkg, 12-place) PLUS any adapter specified in SM-24 tube chart. | | | | | | | | | |
| 16 | 16 | Polysulfone, Flanged Tube | 5250 | 18 x 100 | 03604 (50) | 03269 (50) | Polypropylene, Nonsealing | 00892 (1) | 12 | |
| 16 | 16 | Polypropylene, Flanged Tube | 5250 | 18 x 100 | 03244 (50) | 03269 (50) | Polypropylene, Nonsealing | 00892 (1) | 12 | |
| 16 | 17 | Polycarbonate, Flanged Tube | 5250 | 18 x 100 | 03243 (50) | 03269 (50) | Polypropylene, Nonsealing | 00892 (1) | 12 | |
| 15 | — | Conical, Urinalysis Tube | — | 17 x 120 | — | — | — | 00892 (1) | 12 | |
| 15 | — | Conical, Tissue Culture Tube | — | 17 x 120 | — | — | — | 00892 (1) | 12 | |
| — | — | Chlamydia Vial, Flat Bottomed | — | 15 x 45 | — | — | — | 00726 (1) 00725 (1) | 18 18 | Top Level Adapter 00726 must be used with Bottom Level Adapter 00725. 00725 can be used alone. |
| 15 | — | Vacutainer® Tube | — | 16 x 127 | — | — | — | 00385 (1) | 18 | |
| — | — | BioVial® OmniVial® | — | — | — | — | — | 00725 (1) | 18 | |
| 12 | 12 | Polysulfone, Flanged Tube | 5250 | 16 x 100 | 03602 (50) | 03266 (50) | Polypropylene, Nonsealing | 00385 (1) | 18 | |
| 12 | 12 | Polypropylene, Flanged Tube | 5250 | 16 x 100 | 03116 (50) | 03266 (50) | Polypropylene, Nonsealing | 00385 (1) | 18 | |
| 12 | 12 | Polycarbonate, Flanged Tube | 5250 | 16 x 100 | 03115 (50) | 03266 (50) | Polypropylene, Nonsealing | 00385 (1) | 18 | |

(continued on next page)

TUBES, BOTTLES FOR HG-4L (cont.)

| CAPACITY (ml) | | DESCRIPTION | MAX SPEED (rpm) | DIMENSIONS (mm) | TUBE CAT. NO. (Qty/Pkg) | COVER ASSEMBLY | | ADAPTER REQ'D | | NOTES |
|---------------|------|--|-----------------|-----------------|-------------------------|--------------------|---------------------------|------------------------|---------------|--|
| NOMINAL | FILL | | | | | CAT. NO. (Qty/Pkg) | DESCRIPTION | CAT. NO. (Qty/Pkg) | NO. OF PLACES | |
| 10 | — | Vacutainer® Tube | — | 16 x 100 | — | — | — | 00385 (1) | 18 | |
| 10 | 11 | Polyallomer, Thin Walled, Flanged Tube | 5250 | 13 x 100 | 03124 (50) | — | — | 00384 (1) | 22 | |
| 10 | — | Glass Tube | — | 13 x 100 | — | — | — | 00384 (1) | 22 | |
| 10 | 10 | Pyrex® Tube | 5000 | 13 x 100 | 03118 (50) | — | — | 00384 (1) | 22 | |
| 7 | — | Vacutainer® Tube | — | 13 x 100 | — | — | — | 00384 (1) | 22 | |
| 7 | 7 | Polysulfone, Flanged Tube | 5250 | 13 x 100 | 03601 (50) | 03265 (50) | Polypropylene, Nonsealing | 00384 (1) | 22 | |
| 7 | 7 | Polypropylene, Flanged Tube | 5250 | 13 x 100 | 03121 (50) | 03265 (50) | Polypropylene, Nonsealing | 00384 (1) | 22 | |
| 7 | 7 | Polycarbonate, Flanged Tube | 5250 | 13 x 100 | 03120 (50) | 03265 (50) | Polypropylene, Nonsealing | 00384 (1) | 22 | |
| 5 | 6 | Pyrex® Tube | 5000 | 12 x 75 | 03102 (50) | — | — | 00857 (1) 00858 (1) | 33 33 | Top Level Adapt. 00858 must be used with Bottom Level Adapt. 00857. 00857 may be used alone. |
| 5 | — | Standard Glass Tube | — | 12 x 75 | — | — | — | 00857 (1) 00858 (1) | 33 33 | Top Level Adapt. 00858 must be used with Bottom Level Adapt. 00857. 00857 may be used alone. |
| 5 | — | Vacutainer® Tube | — | 10.25 x 103 | — | — | — | 00857 (1) | 33 | |
| 4 | — | Vacutainer® Tube | — | 10.25 x 82 | — | — | — | 00857 (1) | 33 | |
| 4 | — | Standard Glass Tube | — | 10 x 75 | — | — | — | 00857 (1) 00858 (1) | 33 33 | Top Level Adapt. 00858 must be used with Bottom Level Adapt. 00857. 00857 may be used alone. |
| 4 | 4 | Polypropylene, Flanged Tube | 5250 | 11 x 75 | 03105 (50) | 03264 (50) | Polypropylene, Nonsealing | 00857 (1) 00858 (1) | 33 33 | Top Level Adapt. 00858 must be used with Bottom Level Adapt. 00857. 00857 may be used alone. |
| 4 | 4 | Polycarbonate, Flanged Tube | 5250 | 11 x 75 | 03104 (50) | 03264 (50) | Polypropylene, Nonsealing | 00857 (1) 00858 (1) | 33 33 | Top Level Adapt. 00858 must be used with Bottom Level Adapt. 00857. 00857 may be used alone. |
| 3 | 3 | Pyrex® Tube | 5000 | 10 x 75 | 03100 (50) | — | — | 00857 (1) 00858 (1) | 33 33 | Top Level Adapt. 00858 must be used with Bottom Level Adapt. 00857. 00857 may be used alone. |
| 1.5 | — | Microtube | — | — | — | — | — | 04871 (1) 00857 (1) | — 33 | Adapter plate 04871 must be used together with adapter 00857. |

TO ORDER, CALL 1-800-551-2121